


Optics Lab Data Export Guide

ERAU Virtual Hub

PROCEDURE


STEP 1

Launch the Virtual Hub application by double-clicking on the desktop shortcut.

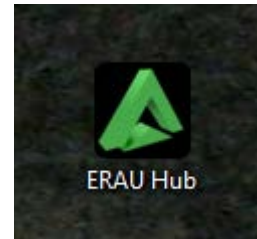


STEP 2

Enter the Optics Lab by pressing E on the door.

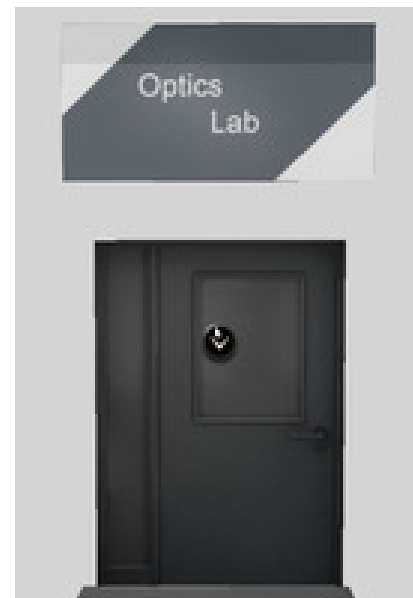

STEP 3

Set the switch to Daylight.

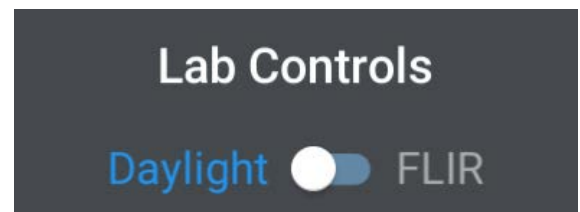

10/24/17



DOC ID



Academic Technology




1.866.554.8406

Page 1

EMBRY-RIDDLE
Aeronautical University.
WORLDWIDE


STEP 4

Set the desired information for the experiment using the sliders.

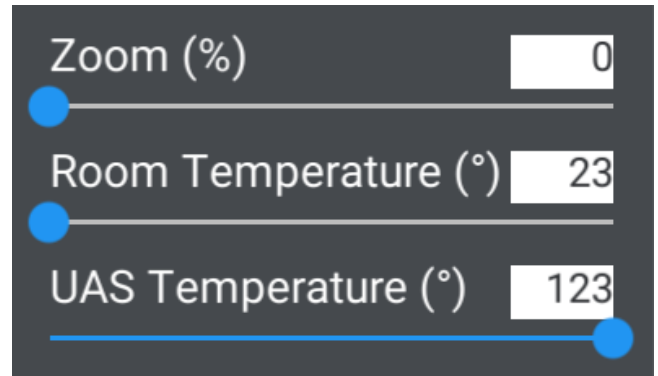


STEP 5

After the temperature information is set move the lab control switch to FLIR.

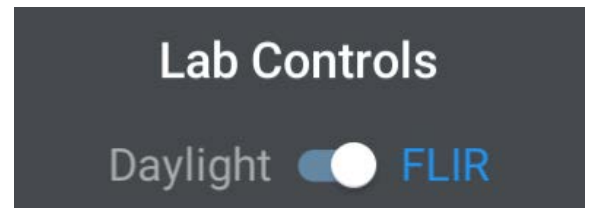

STEP 6

After the experiment, click the “Take Photo” button. The photo will then appear on the photos page. You can access the page using the camera icon on the top left of the screen.

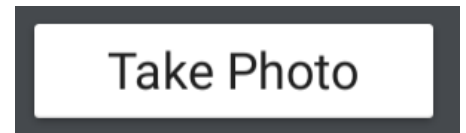



STEP 7

You will select the photos you want then click export. They will appear in a word document with more information about the photo.

Evidence Report

Investigator: User

Photo Name: Photo_2017_08_30_14_52_5
Date Taken: 2017 08 30 14 52 5
Notes:

